

Product Instruction

Leibovitz's L-15 Medium

Product Type: Leibovitz's L-15

Product Description

Leibovitz's L-15 medium contains amino acids, vitamins, inorganic salts, and other nutrients, using phosphate and free basic amino acid buffer system instead of sodium bicarbonate, which is suitable for the primary tissue block cell culture of HEP-2 monkey kidney cells, embryonic and adult tissues. BioEngine has passed the ISO13485 quality management system certification, and the product quality is stable and reliable.

Product Formula

The complete formula can be downloaded from the company's website.

Product Ingredient

The medium contains carbohydrates, amino acids, vitamins, metal ions, and other nutritional components.

This product does not contain hydrolysates, cytokines, antibiotics, HEPES, and nucleosides. This product does not contain raw materials of animal origin.

Product Preservation

- > Store in a dark environment at $2-8^{\circ}$ C.
- This product is vulnerable to water damage. Please use it immediately after opening. If it needs to be stored, please seal it by heat sealing and sealing clips, avoiding dampness and being ineffective.
- > Do not recommend using when the product is beyond the expiration date.

Instructions for the preparation of Leibovitz's L-15 medium

Prepare Leibovitz's L-15 medium, as Table 1 shows.

onent	Concentration
eibovitz's L-15	14.08g/L
um	
um	

Table 1 Preparation of Leibovitz's L-15 medium

Shanghai BioEngine Sci-Tech Co., Ltd.

5F, Building 1, Lane 720, Cailun Road, Pudong New Area, Shanghai Tel: 021-68582660 www.bio-engine.com.cn



- (1) Weigh <u>90%</u> water <u>of</u> the final medium <u>preparation volume</u> into the medium preparation container. When preparing, ultrapure water or water for injection and above standard water should be used, and the water temperature should be controlled at 20-30°C. Turn on the mixing system of the medium preparation container, stir thoroughly, and avoid the generation of air bubbles during stirring.
- (2) Accurately weigh the corresponding mass of Leibovitz's L-15 medium dry powder according to the concentration ratio of 14.08g/L, add it to the preparation container of (1), and stir thoroughly for 15-30 minutes until it becomes clear.
- (3) Use 5mol/L sodium hydroxide solution or 1mol/L hydrochloric acid solution to adjust the pH value of the medium to 7.60, and use preparation water to <u>quantify to 100% of the preparation volume</u>.
- (4) It is recommended to use a pulse pump or compressed air (3-15 psi) to sterile filter the Leibovitz's L-15 medium solution through a sterile filter membrane with 0.22 μm pore size.
- (5) The prepared medium liquid should be stored in a dark environment at 2-8°C.
- (6) Reference parameters of the final product

Indicator	Reference Standard
Product pH value	7.30-7.90
Osmolality	290-340mOsm/Kg

Notes:

- (1) The above units of "g/L" are volume concentration (solute mass/solution volume).
- (2) The above preparation parameters (such as stirring time, etc.) are for small-scale preparation in research and development. When in mass production, please set appropriate preparation parameters according to the stirring capacity of the preparation vessel.