



Proli-S001S PK15 Serum-Free Media

A serum-free medium especially developed for PK15 cell culture and efficient amplification of circovirus

Proli-S001S is an innovative serum-free medium developed independently by Shanghai BioEngine Sci-Tech Co., Ltd., which is serum-free, protein-free and animal-derived component-free. It is suitable for PK15 suspension cell culture. It supports efficient acclimation of adherent PK15 cells and high density growth of suspension cells, and provides a high quality environment for efficient amplification of porcine circovirus. *Proli-S001S* is your ideal choice for large-scale industrial suspension culture of PK15 cells: the suspension culture system provides you with a significant increase in production efficiency, while the serum-free culture system provides you with maximum process stability. It has been industrially applied in the production of porcine circovirus vaccine project.

Features

- Serum-free
- S Animal-derived component-free
- Protein-free
- Support rapid serum-free suspension adaptation of adherent PK15 cells
- Supporting high-density growth of PK15 cells
- Supporting PK15 cells for efficient production of circovirus

Advantages

- Solution Full traceability by EU-certified ISO13485:2016 Quality anagement System;
- Second Se
- Optional powder media for use in large-scale manufacturing with easy preparation procedures;
- Powder media capable of a single batch size of 100,000 L.

*CPK: Process Capability Index; a CPK>1.33 indicates good process control and small inter-batch difference in products

Ordering Information

Product Name	Cat. No.	Form	Size	Package
Proli-S001 PK15 Serum-free Medium	EXP0101001	Liquid	1L	Bottle
Proli-S001S PK15 Serum-free Medium	EXP0111103	Powder	10L	Bag
	EXP0111102	Powder	100L	Bag
	EXP0111101	Powder	200L	Bag



Proli Series PK15 Serum-Free Medium


Performance

Scell adaptation

Adherent PK15 cells were directly acclimated in Proli-S001S and got stabilized after 10 generations, and the specific growth rate gradually increased with the number of passages, with a doubling time of 22-25h after stabilization.

Sell growth

Proli-S001S supports high density growth of suspension PK15 cells up to 107 cells/ml.

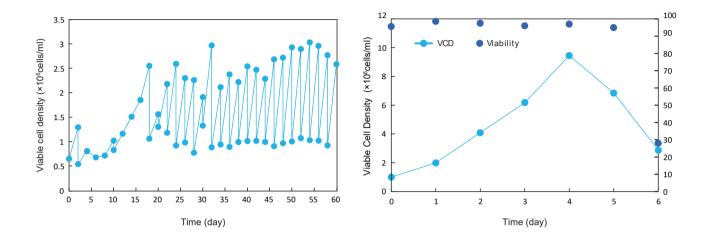


© Virus production

Add.:


0

Using Proli-S001S for culture and suspension of PK15, either for (a) fluorescence quantitative PCR or (b) TCID50 assay, showed substantially higher titers of virus production compared to the original adherent process.

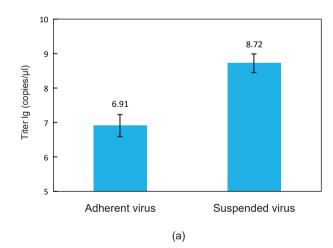

Floor 5, Building 1, Lane 720, Cailun Road, Zhangjiang

Hi-Tech Park, Pudong, Shanghai, PRC

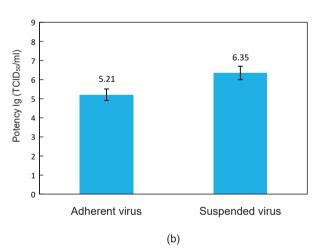

30 years of ingenuity on creating a novel drive for cell culture



Tel.: (86)21-68582660

Web .: Web.: www.bio-engine.com.cn

 \mathbf{M}

